

Technical Article

Chocolate Challenge! See How Spectroscopy Can Give You Answers.



Stuti Ghiya

Roses are red,

Violets are blue,

We used a TI DLP® NIRscan™ Nano EVM,

To pick the right chocolate for you.

Love may be fleeting, but at least we'll always have chocolates.

While there's nothing sweeter for your sweetheart than a box of chocolates on Valentine's Day, choosing the option with just the right balance of cocoa and sugar content for your taste buds can be a lovelorn experience.



Figure 1. The DLP NIRscan Nano EVM in action.

At TI DLP Products, we decided to address this candy conundrum with our [DLP NIRscan Nano evaluation module \(EVM\)](#). We scanned unidentified samples of chocolate and compared it to a database of known milk chocolates and dark chocolates of approximately 70 percent cocoa and 85 percent cocoa.

Based on the known chocolates in the database, here's what our testing found:

- 70 percent dark chocolate = approximately 27.5 percent sugar
- 85 percent dark chocolate = approximately 12.5 percent sugar
- Milk chocolate = approximately 52.4 percent sugar

The DLP NIRscan Nano EVM is optimized for a 900 to 1,700 nanometer wavelength range and features a 6,000:1 signal-to-noise ratio. The results shown below were taken directly from the EVM's included graphical user interface (GUI). We observed the absorbance levels of all three varieties of chocolate across this wavelength range. When compared to the absorbance spectrums in the database, we could identify which chocolates were 70 percent cocoa, 85 percent cocoa, or milk chocolate.


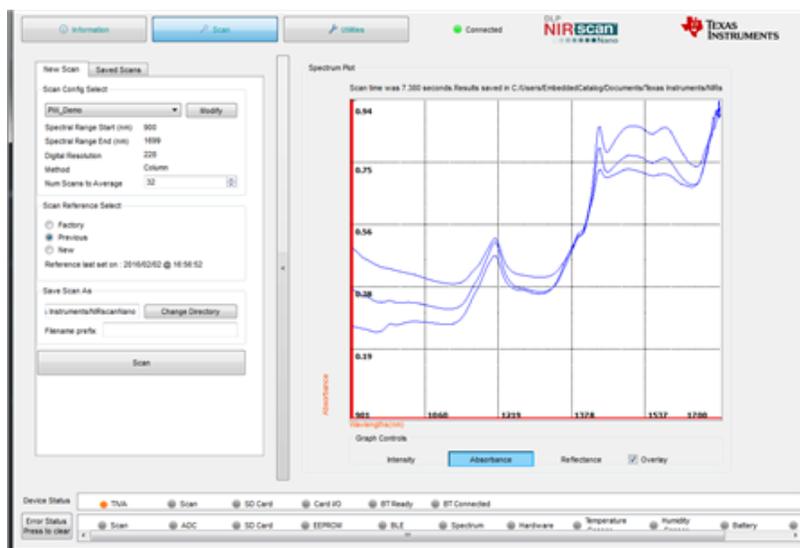


Figure 2. This screen shot of the GUI plots the absorbance results of the three types of chocolate we tested.

So Much More than a Box of Chocolates

Of course, the NIRscan Nano EVM can help determine more than just sugar content.

The EVM is an evaluation tool to help developers quickly design spectrometers and chemical analyzers using DLP chips. It uses our DLP technology in conjunction with a single element InGaAs detector to deliver high performance measurements in a portable form factor that is more affordable than expensive InGaAs array detectors or fragile rotating grating architectures. The wide NIR-window transmittance also makes it useful in a variety of industries such as agriculture, petrochemicals, and health care.

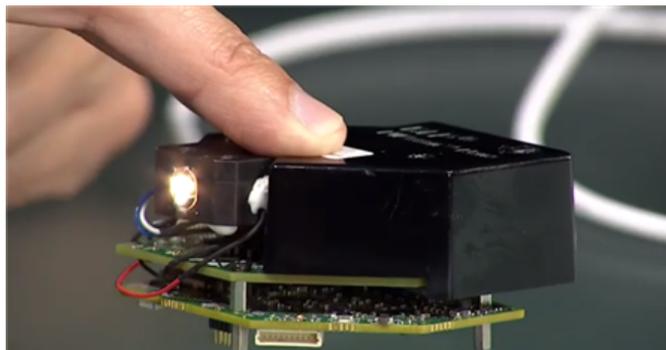

[DLP NIRscan Nano EVM: Getting Started Out of the Box](#)

The design can be powered via USB or batteries and is packed with TI components, including the [DLP2010NIR](#) digital micromirror device (DMD), [DLPC150](#) digital controller, and [DLPA2005](#) power management and LED/lamp driver. The DLP2010NIR chip is our smallest, most efficient DLP chip, designed for handheld near-infrared sensing application areas.

Now you can identify the right chocolate for your taste buds from a box of chocolates and know just how much sugar these three types of chocolate contain. How sweet is that?

IMPORTANT NOTICE AND DISCLAIMER

TI PROVIDES TECHNICAL AND RELIABILITY DATA (INCLUDING DATA SHEETS), DESIGN RESOURCES (INCLUDING REFERENCE DESIGNS), APPLICATION OR OTHER DESIGN ADVICE, WEB TOOLS, SAFETY INFORMATION, AND OTHER RESOURCES "AS IS" AND WITH ALL FAULTS, AND DISCLAIMS ALL WARRANTIES, EXPRESS AND IMPLIED, INCLUDING WITHOUT LIMITATION ANY IMPLIED WARRANTIES OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE OR NON-INFRINGEMENT OF THIRD PARTY INTELLECTUAL PROPERTY RIGHTS.

These resources are intended for skilled developers designing with TI products. You are solely responsible for (1) selecting the appropriate TI products for your application, (2) designing, validating and testing your application, and (3) ensuring your application meets applicable standards, and any other safety, security, regulatory or other requirements.

These resources are subject to change without notice. TI grants you permission to use these resources only for development of an application that uses the TI products described in the resource. Other reproduction and display of these resources is prohibited. No license is granted to any other TI intellectual property right or to any third party intellectual property right. TI disclaims responsibility for, and you will fully indemnify TI and its representatives against, any claims, damages, costs, losses, and liabilities arising out of your use of these resources.

TI's products are provided subject to [TI's Terms of Sale](#) or other applicable terms available either on [ti.com](https://www.ti.com) or provided in conjunction with such TI products. TI's provision of these resources does not expand or otherwise alter TI's applicable warranties or warranty disclaimers for TI products.

TI objects to and rejects any additional or different terms you may have proposed.

Mailing Address: Texas Instruments, Post Office Box 655303, Dallas, Texas 75265
Copyright © 2023, Texas Instruments Incorporated